07553-77 with pump head 77201-60

OPERATING MANUAL:
PUMP DRIVES

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A-1299-0637
Edition 03

SAFETY PRECAUTIONS

DANGER: High voltages exist and are accessible in the Drive/Controller.

WARNING: Disconnect the AC power input line cord before connecting the drive motor cable.

WARNING: No user serviceable parts are inside of this controller. Refer servicing to your dealer.

WARNING: PRODUCT USE LIMITATION

This product is not designed for, nor intended for use in patient connected applications; including, but not limited to medical and dental use and accordingly, has not been submitted for FDA approval.

Variety of Pump Heads Accepted

Mount up to 2 (600 rpm) or 4 (100 rpm) MASTERFLEX® L/S® Pump Heads and all MASTERFLEX-compatible pump heads.

Setup and Drive Operation

1. Mount pump head and load tubing.
(See pump head manual.)
2. Connect the drive motor cable to the receptacle on the rear side of the controller.
3. Connect the AC power input line cord to the AC receptacle.
4. Selecting motor direction turns pump on.
5. Adjust flow rate with speed control potentiometer.

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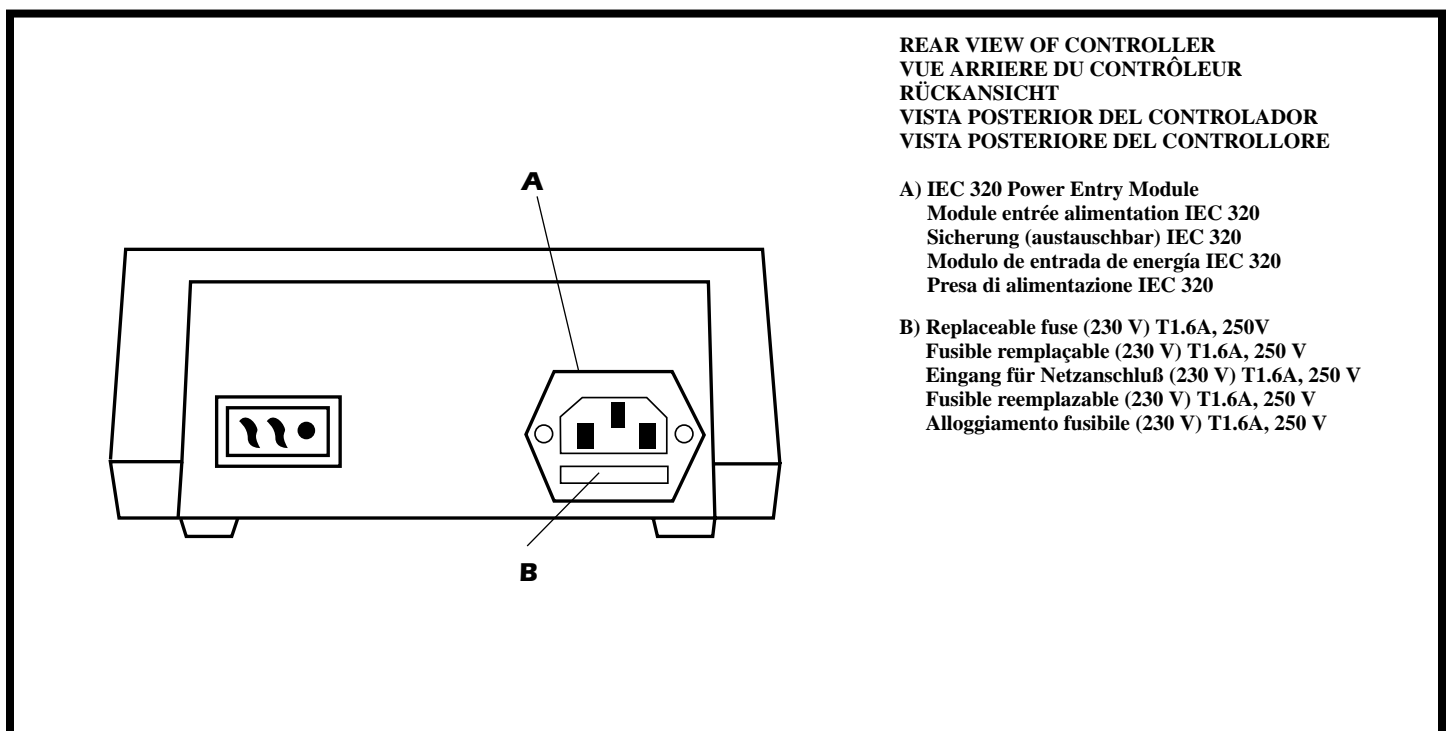
MOTOR BRUSH CHECK/REPLACEMENT

NOTE: Brushes should be checked every 6 months or 3000 operating hours.

1. Place the FWD-OFF-REV switch in the OFF position.
2. Disconnect the AC line cord from the AC receptacle.
3. Disconnect the drive motor cable from the receptacle on the rear of the controller.
4. Carefully unscrew each brush holder on opposite sides of the motor. Withdraw the brush, and examine it for wear.

NOTE: Replace both brushes, if either brush is less than 0.300 in long from base to point.

5. Screw brushes into brush holder on each side of motor.
6. Connect motor to rear panel connector on controller.
7. Connect AC line cord to AC receptacle.

Fuse Replacement

1. On 230V units, place the FWD-OFF-REV switch in the OFF position.
2. Disconnect the AC line cord from the rear of the unit.
3. Remove and check the fuse located on the rear panel below the AC power connector and replace if defective.
4. Reconnect AC line cord.

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Name of Apparatus: MASTERFLEX® Analog Modular Benchtop L/S® Drive

Model Numbers: 7553-77, 7553-78*, 7553-79*, 7553-87, 7553-89*, 77913-25

Description of Apparatus: Variable Speed, Peristaltic Pump Motor Drive. Used with peristaltic pump head and tubing to pump fluids.

7553-77 – Modular System, 600 rpm

7553-79 – Modular Motor, 600 rpm

7553-87 – Modular System, 100 rpm

7553-89 – Modular Motor, 100 rpm

7553-78 – Modular Controller

77913-25 – Bundled 7553-77

Barnant Company declares that the above models are in conformity to the following harmonized standards and directives:

Applicable Directives	Applicable Specifications	Manufacturer's Report Number
73/23/EEC 93/68/EEC	EN61010-1/A2: 1995	TR9538
89/336/EEC 92/31/EEC 93/68/EEC	EN61326-1/A1: 1998	TR9539

The last two digits of the year in which the current configuration of the above models was assessed per the Low Voltage Directive is: 00.

*Evaluated for EMC as part of a system.

Manufacturer: Barnant Company Division
Cole-Parmer Instrument Company
28W092 Commercial Avenue
Barrington, IL 60010-2392
USA
Tel.: 847-381-7050

Manufacturer's Signature:

James W. Doll
Vice President, Engineering

5 October, 2000

Date

WARRANTY

Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is noted on the Warranty Card.

PRODUCT RETURN

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

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Vernon Hills, Illinois U.S.A. 60061-1844
Barrington, Illinois U.S.A. 60010-2392