7592-90 Wall Mount Controller with 7592-40 Pump Drive
and 7592-82 Benchtop Controller

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Name of Apparatus: MASTERFLEX® Digital Modular I/P® Pump Drive

Model Numbers: 7592-27, -35, -45*, -85* and -95*, 77970-27 and -37

Description of Apparatus: Digital-controlled, Variable Speed, Peristaltic Pump and Drive. Used with peristaltic tubing to pump fluids.

- 7592-27 – Modular Benchtop System
- 7592-35 – Modular NEMA System
- 7592-85 – Modular Benchtop Controller
- 7592-95 – Modular NEMA Controller
- 7592-45 – Modular Motor
- 77970-27 – Bundled 7592-27
- 77970-37 – Bundled 7592-35

Barnant Company declares that the above models are in conformity to the following harmonized standards and directives:

Applicable Directives	Applicable Specifications	Manufacturer's Report Number
73/23/EEC 93/68/EEC	EN61010-1/A2: 1995	TR9864
89/336/EEC 92/31/EEC 93/68/EEC	EN61326-1/A1: 1998	TR9865

The last two digits of the year in which the current configuration of the above models was assessed per the Low Voltage Directive is: 00.

*Evaluated for EMC as part of a system.

Manufacturer: Barnant Company Division
 Cole-Parmer Instrument Company
 28W092 Commercial Avenue
 Barrington, IL 60010-2392
 USA
 Tel.: 847-381-7050

Manufacturer's Signature:

James W. Doll
 James W. Doll
 Vice President, Engineering

16 August, 2000
 Date

PUMP DRIVES—MODEL NOS.

TYPE	SYSTEM	PUMP DRIVES	CONTROLLER
Benchtop (115V)	7592-20	7592-40	7592-82
Benchtop (230V)	7592-27	7592-45	7592-85
NEMA (115V)	7592-30	7592-40	7592-90
NEMA (230V)	7592-35	7592-45	7592-95

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TYPE	SYSTÈME	ENTRAÎNEMENTS DE POMPES	CONTRÔLEUR
Banc (115 V)	7592-20	7592-40	7592-82
Banc (230 V)	7592-27	7592-45	7592-85
NEMA (115 V)	7592-30	7592-40	7592-90
NEMA (230 V)	7592-35	7592-45	7592-95

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SAFETY PRECAUTIONS

DANGER: High voltages exist and are accessible in the Modular Drive. Use extreme caution when servicing internal components.

WARNING: Tubing breakage may result in fluid being sprayed from pump. Use appropriate measures to protect operator and equipment.

WARNING: Turn drive off before removing or installing tubing. Fingers or loose clothing could get caught in drive mechanism.

WARNING: PRODUCT USE LIMITATION

This product is not designed for, nor intended for use in patient connected applications; including, but not limited to, medical and dental use, and accordingly has not been submitted for FDA approval.

Introduction

The Modular Drive controls the speed of MASTERFLEX® I/P® (Industrial/Process) pumps to provide flow rates from 0.12 to 13 L/min.

Mount up to 2 MASTERFLEX® I/P® pump heads and all MASTERFLEX-compatible pump heads.

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Control/Display Functions/ Mounting Dimensions

Press buttons to activate function.

Use up/down (▲, ▼) arrows to correct/change a flashing display.

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A) DOWN ARROW (DECREMENT)—Decrease value of a flashing display.

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B) UP ARROW (INCREMENT)—Increase value of a flashing display.

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C) DISPENSE/COPY—Select dispensed volume, number of copies of a dispensed volume, or the dispense interval.

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D) FLOW CONTROL—Set flow rate for selected tubing size. To change flow rate, press ▲ or ▼ arrows. (If pump is running, its speed will change with new settings.)

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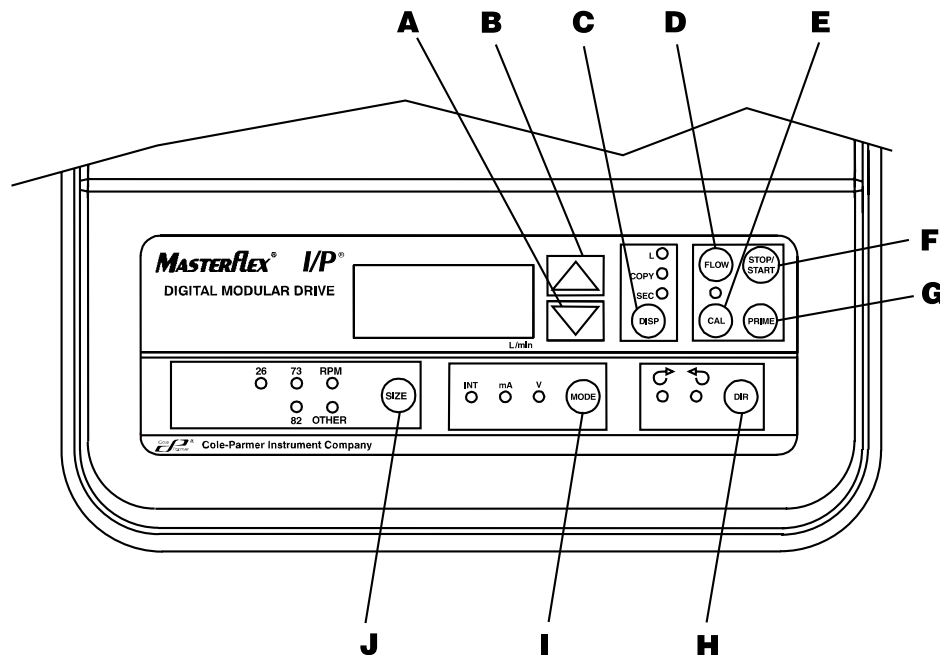
E) CAL CONTROL—Refine built-in calibration, using a measured volume.

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CONTROL/DISPLAY FUNCTIONS/MOUNTING DIMENSIONS

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G) PRIME—Run pump at full speed to fill or clear lines.

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H) DIRECTION—To change motor direction.

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I) MODE SELECT—INT for internal control; mA for remote current control; V for remote voltage control.

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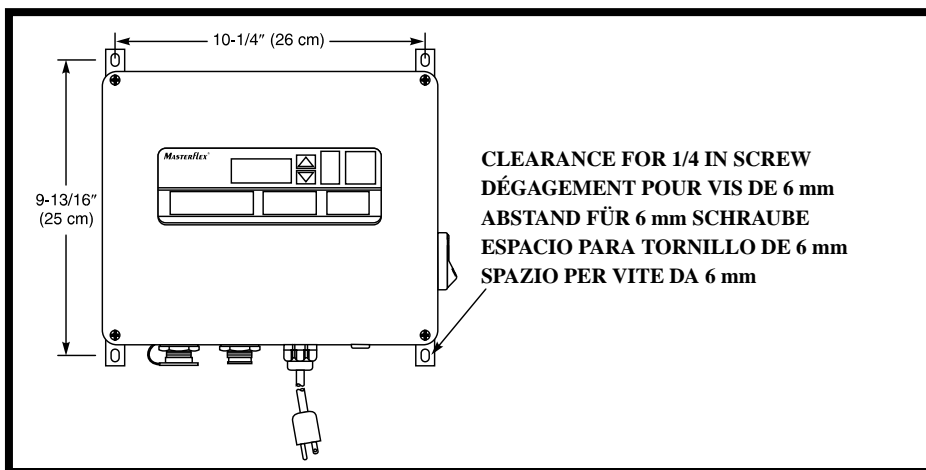
Mounting Dimensions for Model Numbers 7592-30, -35

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Setup and Drive Operation

1. Connect Motor Cable plug to mating receptacle on the Controller.
2. Connect power cord to Controller and grounded power line outlet.
3. Mount pump head and load tubing. (See pump head manual.)
4. Turn pump on and select TUBING SIZE.

NOTE: If CAL LED is lit, that tubing size has been previously field calibrated. If LED is not lit, the drive is operating with the built-in factory calibration. To clear a field calibration, press and hold the CAL switch until the CAL light goes out. This will take about 3 seconds. To recalibrate for better accuracy, see Calibration section.

5. MODE selection (INT, mA, V).
6. Select MOTOR DIRECTION (CW or CCW).
7. PRIME and CALibrate the pump (if required).
8. Press FLOW button and watch display to set the flow rate with UP/DOWN keys.

9. Press STOP/START button to begin pumping.

NOTE: Pump will restart automatically after a brownout or powerout condition.

Calibration

Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

1. Select correct tubing size and flow rate.
2. Press CAL, calibration volume appears.
3. Press STOP/START, the pump will use its stored memory to dispense the specified calibration sample quantity. The pump will stop automatically.
4. Weigh/measure the sample.
5. Use UP/DOWN arrow keys to correct the flashing display.

NOTE: If the adjusted calibration is too great, "Err" will appear in the display. If this occurs, press the CAL control and repeat the calibration procedure. The microprocessor will retain one special calibration value per tubing size, even when power is turned off. The next calibration will replace the existing value.

6. Press SIZE to exit the calibration cycle.

Maximum Flowrate (OTHER Tubing)

1. To set the maximum flowrate for non-standard pumps or tubing sizes OTHER, press CAL, then FLOW. The maximum flowrate will then flash on the display.
2. Use UP/DOWN arrow keys to set desired flowrate.
3. Press SIZE to exit.

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SETUP AND DRIVE OPERATION CALIBRATION

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DISPense/copy

A first press of the DISP key results in the last entered dispense volume being displayed. The "L" annunciator will illuminate and flash. The INC/DEC keys are used to change the dispense volume, if desired. The STOP/START key then initiates delivery of the set volume. The amount remaining to be dispensed will be displayed during countdown. The dispense function is exited by pressing any key except Increment, Decrement, DISP, or STOP/START.

A second press of the DISP key causes the COPY annunciator to illuminate and flash. The STOP/START key is then used to deliver the desired volume without the need to know the volume in specific units. A third press of the DISP key enters the volume dispensed. The COPY annunciator stops flashing. The STOP/START key is then used to initiate delivery of the copied volume. The number of copies dispensed will be displayed after each dispense. The maximum number of copies is 9999. The STOP/START key is used to pause the copy dispense during dispensing; copy dispense can then be continued using the STOP/START key.

A fourth press of the DISP key results in the last entered dispense interval being displayed. The SEC annunciator will illuminate and flash. The INC/DEC keys are used to change the dispense time, if desired, from 1 to 9999 seconds. The STOP/START key then initiates delivery for the set time interval. The remaining time will be displayed during countdown. Pressing the DISP key a fifth time exits this mode.

Keypad Lockout Enable/Disable

Press and hold FLOW. After five (5) seconds, display will change to all dashes. Then, while holding FLOW, press PRIME five (5) times.

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REMOTE CONTROL

Selectable input (0–20 mA, 4–20 mA, 0–10V DC)
 ±0.5% linearity control
 2300V isolation potential
 START/STOP; CW/CCW; PRIME via contact closure

Remote Control Setup

1. Place the power switch in the off position.

CAUTION: Power must be turned off before connecting the external remote control cable to prevent damage to the drive.

2. Connect the cable from the external remote control to the DB15-pin receptacle on the rear panel. On washdown units, connect to the mating receptacle on the bottom panel.

3. Select type of remote control input and output required as follows:

- a). Press and hold the MODE control while turning the power switch to the on position. After two seconds, release the MODE control. The initial display will show: "inP". After two seconds the display will show either 0–20 or 4–20.

NOTE: Press the up (increment) or down (decrement) arrows to select between 4–20 and 0–20 for current loop control.

- b). Press the MODE control again. The initial display will show: "out". After two seconds the display will show either 0–20 or 4–20.

NOTE: Press the up (increment) or down (decrement) arrows to select between 4–20 and 0–20 for current loop output.

4. Press the MODE control to select mode of operation. The LED's indicate the selected mode. Select either mA or V.

NOTE: If only remote STOP/START, PRIME and/or CW/CCW is to be used, the MODE control can be set to any of the three positions.

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REMOTE CONTROL COMMANDE À DISTANCE EXTERNE STEUERUNG CONTROL REMOTO TELECOMANDO

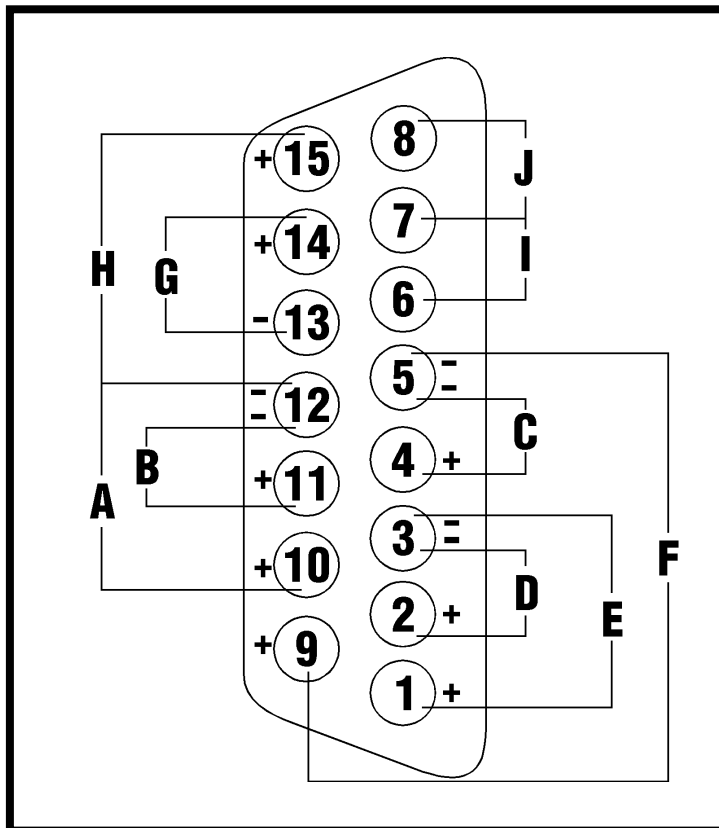
Connector pin configuration/with wiring scheme

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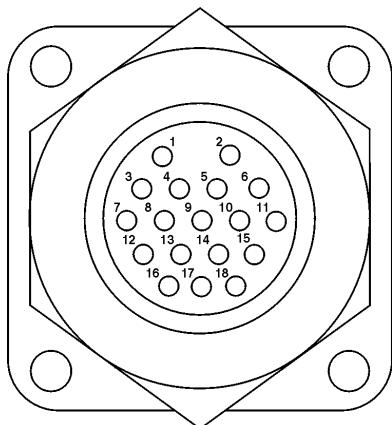
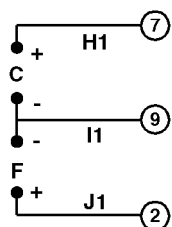
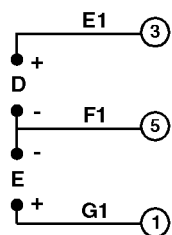
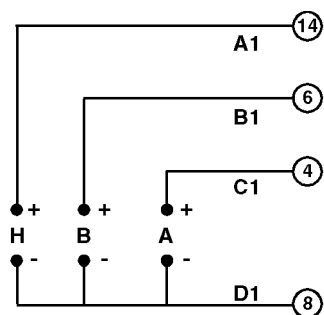
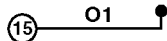
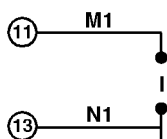
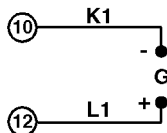
NOTE: Colors are those of Remote Cable, Cat. number MN-77300-32.

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TROUBLESHOOTING AND MAINTENANCE

DANGER: Modular Drive Systems are not customer serviceable, except for replacement of FUSES and MOTOR BRUSHES.

WARNING: Unplug Controller prior to checking or replacing FUSES and/or MOTOR BRUSHES.

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A) T6.3A FUSE (115V AC); T3.15A FUSE (230V AC)
CAUTION: Do not substitute.

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D) MOTOR RECEPTACLE

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RECEPTÁCULO DE MOTOR

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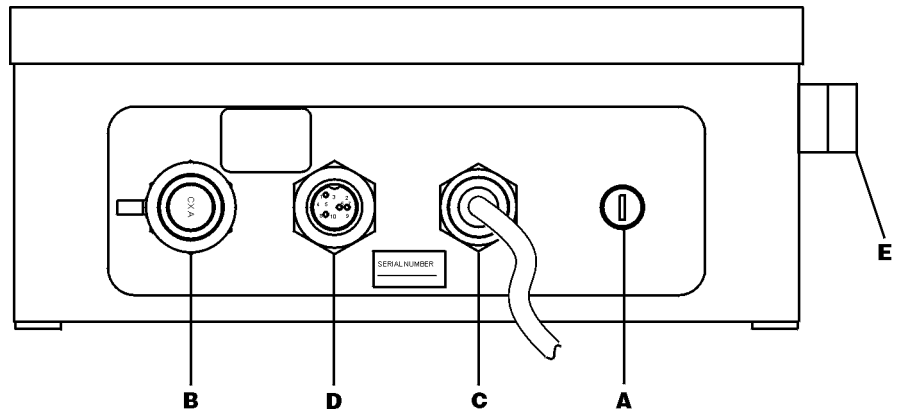
E) POWER SWITCH — ALL SETTINGS ARE RETAINED IN MEMORY

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7592-90, -95

MOTOR BRUSH CHECK/REPLACEMENT

Brushes should be checked every 6 months or 2000 operating hours or if erratic operation occurs.

Erratic operation may occur immediately after brush replacement. Allow motor to run up to an hour to allow brushes to seat.

DANGER: High voltages exist and are accessible in the Modular Drive. Other than replacement of FUSES and/or MOTOR BRUSHES, refer servicing to qualified personnel.

To replace the brushes, remove the 2 rubber caps and unscrew the brushes.

Inspect for cracks and excessive wear.

Replace both brushes, if either brush is less than 7.6 mm (0.300 in) long from base to point.

Contact your dealer if you have service needs.

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Cleaning

Keep the drive enclosure clean with mild detergents. Never immerse nor use excessive fluid.

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SPECIFICATIONS	
Output:	
Speed:	20 to 650 r/min
Torque output, Maximum:	300 oz-in (22 kg•cm)
Speed regulation:	
Line	±0.25% F.S.
Load	±0.25% F.S.
Drift	±0.25% F.S.
Display:	Four-digit, seven segment LED
Remote outputs:	
	Voltage speed output (0–10V DC)
	Current speed output (0–20 mA or 4–20 mA)
	Tach output (TTL, 128 to 7680 Hz)
	Motor running output (N.O. & N.C. contact closure)
Input:	
Supply voltage limits:	
07592-20, -30	90 to 130 Vrms @ 50/60 Hz
07592-27, -35	200 to 260 Vrms @ 50 Hz
Current, max.:	
07592-20, -30	4.4A
07592-27, -35	2.2A
Installation Category:	Installation Category II per IEC 664 (Local level — appliances, portable equipment, etc.)
Remote Inputs:	
	START/STOP, CW/CCW, PRIME (Contact closure)
	Voltage input (0–10V DC)
	Current input (0–20 mA or 4–20 mA)
Construction:	
Dimensions (L × W × H):	
07592-82, -85	9 3/4 in × 9 3/16 in × 5 1/16 in (248 × 233 × 129 mm)
07592-90, -95	9 in × 11 in × 4 1/2 in (229 × 279 × 114 mm)
07592-40, -45	14 1/8 in × 5 9/16 in × 6 in (359 × 141 × 152 mm)
Weight:	
07592-82, -85	4.0 lbs (1.8 kg)
07592-90, -95	9.4 lbs (4.3 kg)
07592-40, -45	16.3 lbs (7.4 kg)
Enclosure Rating:	
07592-82, -85	IP 22 per IEC 529
07592-90, -95	IP 56 (NEMA 4) per IEC 529
07592-40, -45	IP 34 (NEMA 4) per IEC 529
Environment:	
Temperature, Operating:	0° to 40°C (32° to 104°F)
Temperature, Storage:	–45° to 65°C (–49° to 149°F)
Humidity (non-condensing):	10% to 90%
Altitude:	Less than 2000 m
Pollution Degree:	
07592-20, -27	Pollution Degree 2 per IEC 664 (Indoor use — lab, office)
07592-30, -35	Pollution Degree 3 per IEC 664 (Sheltered locations)
Chemical Resistance:	Exposed material is painted aluminum, plastic and vinyl
Compliance:	
	115V: UL 508, CSA C22.2, No. 14-M91
	230V (For CE Mark): EN61010-1/A2: 1995 (EU Low Voltage Directive) and EN61326-1/A1: 1998 (EU EMC Directive)

**REPLACEMENT PARTS
AND ACCESSORIES;
CLEANING;
SPECIFICATIONS**

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The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is noted on the Warranty Card.

PRODUCT RETURN

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

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Vernon Hills, Illinois U.S.A. 60061-1844
Barrington, Illinois U.S.A. 60010-2392