OPERATING MANUAL:
PUMP DRIVES

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EU Declaration of Conformity

Name of Apparatus: MASTERFLEX[®] Console Analog L/S[®] Pump Drives

Model Numbers: 7520-47, -57, -67; 77521-47, -57; 77911-27; 77912-07; 77914-07; 77915-07

Description of Apparatus: Variable speed, peristaltic pump motor drive. Used with pump head and peristaltic tubing to pump fluids.

7520-47 (600 rpm), -57 (100 rpm), -67 (300 rpm)

77521-47 (600 rpm w/Remote), -57 (100 rpm w/Remote);

77911-27 & 77915-07 (Bundled 77521-47);

77912-07 (Bundled 7520-67);

77914-07 (Bundled 7520-57);

Barnant Company declares that the above models are in conformity to the following harmonized standards and directives:

Applicable Directives	Applicable Specifications	Manufacturer's Report Number
73/23/EEC 93/68/EEC	EN61010-1/A2:1995	TR0071
89/336/EEC 92/31/EEC 93/68/EEC	EN61326-1/A1:1998	TR0072

The last two digits of the year in which the current configuration of the above models was assessed per the Low Voltage Directive is: 00.

Manufacturer: Barnant Company Division
Cole-Parmer Instrument Company
28W092 Commercial Avenue
Barrington, IL 60010-2392
USA
Tel: 847-381-7050

Manufacturer's Signature:

James W. Doll
Vice President, Engineering

23 June, 2000

Date

SAFETY PRECAUTIONS

DANGER: High voltages exist and are accessible in the Console Drive. Use extreme caution when servicing internal components.

WARNING: PRODUCT USE LIMITATION

These products are not designed for, nor intended for use in patient connected applications; including, but not limited to, medical and dental use, and accordingly have not been submitted for FDA approval.

Variety of Pump Heads Accepted

Mount up to 2 (300 and 600 rpm) or 4 (100 rpm) MASTERFLEX® L/S® Pump Heads and all MASTERFLEX compatible pump heads.

Setup and Drive Operation

1. Mount pump head and load tubing. (See pump head manual.)
2. Set Mode Select switch (LOCAL/REMOTE).
3. Selecting motor direction turns pump on.
4. Adjust flow rate with the 1- or 10-turn potentiometer speed control.

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REMOTE OPERATING PROCEDURE

WARNING: Turn off AC power before connecting the external remote control cable.

Description

Models with remote inputs (77521 series) have a 10-turn speed control pot with turns counting dial, which replaces the single-turn speed pot. The drive also incorporates a rear panel mounted LOCAL/REMOTE switch along with a 9-pin "D" shell connector (see ACCESSORIES, 07595-45) for an external 2–10V DC or 4–20 mA Speed Control input and Remote Start/Stop operation from external contact closures. This connector incorporates the following Feature Controls, (shown below in Figure 1).

Remote Start/Stop Control

The LOCAL/REMOTE selector switch (located on rear panel) selects the speed control signal from either the front panel or the Remote Speed Control. The Remote Start/Stop function is also enabled or disabled through this selector switch. The Remote/Local contact closure input is only active in the external mode of operation, allowing a remote selection of the speed control command from either the linear inputs or the front panel speed pot. An active low enables the front panel speed pot. (Footswitch available, 77595-35 as listed in the ACCESSORIES.)

The Remote Start/Stop operates through either an open contact closure or open collector driver. An active low causes the motor to run.

The table on page 8 summarizes the remote function operation.

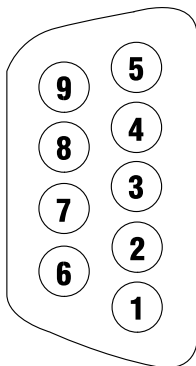
Figure 1
Abb. 1
Figura 1

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
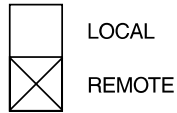
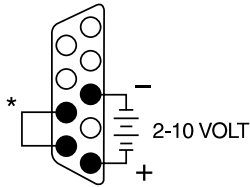
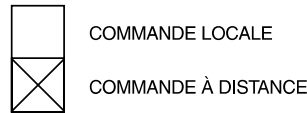
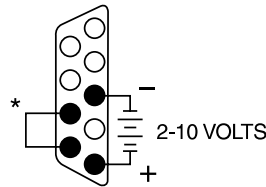
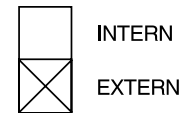
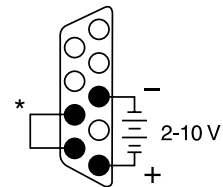
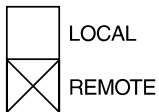
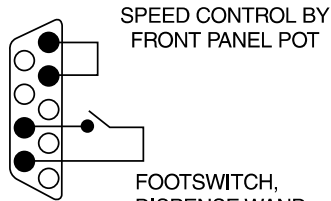
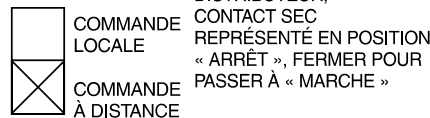
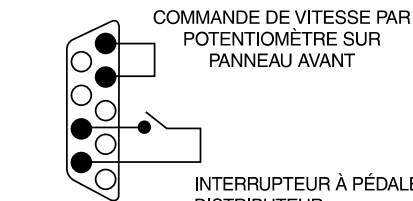
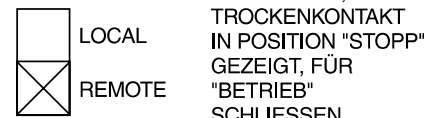
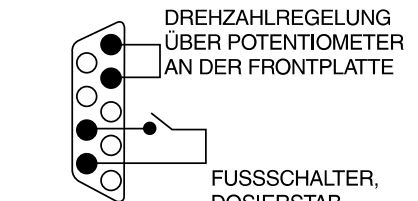
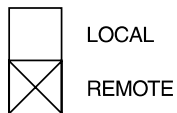
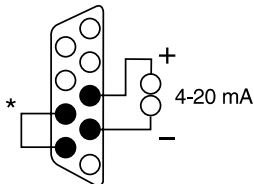
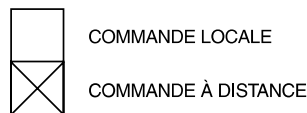
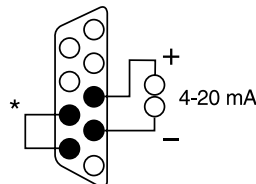
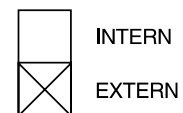
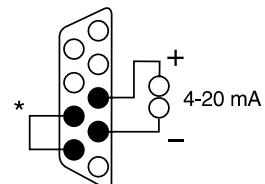
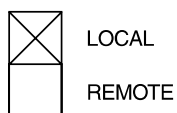
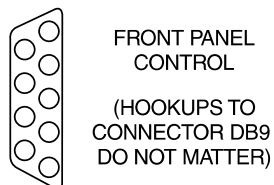
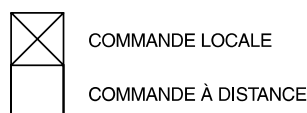
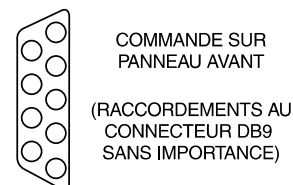
EXEMPLE 4

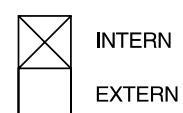
BEISPIEL 4

* JUMPER CAN BE REPLACED WITH SWITCH SHOWN IN EXAMPLE 2

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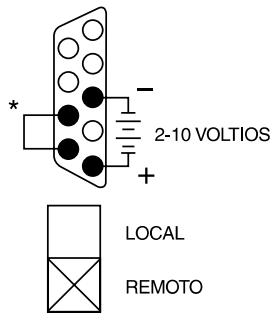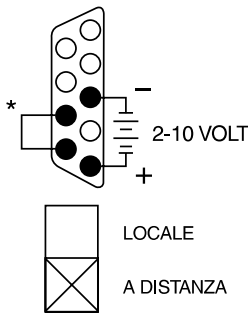
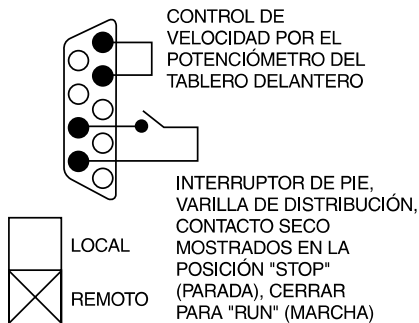
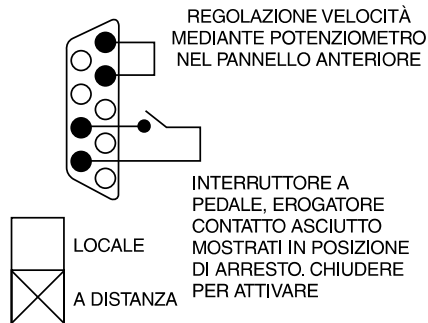
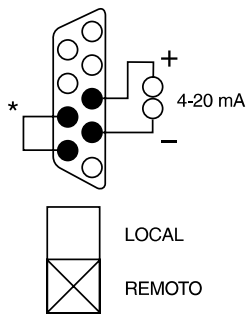
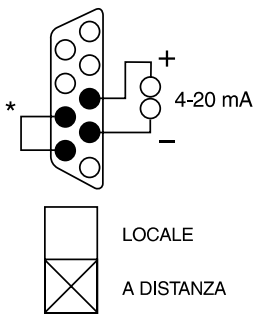
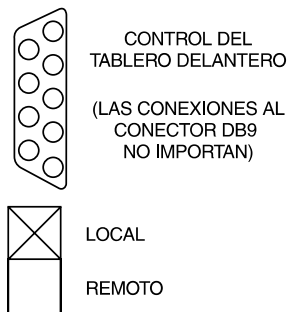
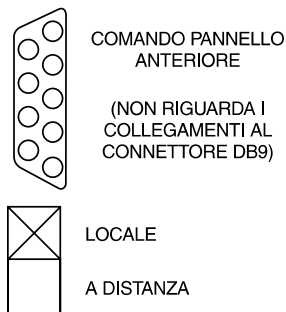
MOTOR BRUSH CHECK/REPLACEMENT

NOTE: Brushes should be checked every 6 months or 3000 operating hours or if erratic operation occurs.

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1. Place the POWER/DIRECTIONAL SWITCH in the off position.
2. Disconnect the AC power input line cord from the AC receptacle.
3. Remove the three screws from each side of the housing and lift off the housing.
4. Carefully unscrew each brush cap. Withdraw the brush, and examine it for wear.

NOTE: Replace both brushes, if either brush is less than 7.6 mm (0.300 in) long from base to point.

5. Insert brushes and install brush cap.
6. Install housing and secure with the three screws on each side.
7. Connect the AC power input line cord to the AC receptacle.

Fuses

NOTE: These drives have an External fuse; T3.15A (115V AC); T1.6A (230V AC). (See REPLACEMENT PARTS / ACCESSORIES Section.)

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MOTOR BRUSH CHECK/REPLACEMENT FUSES CLEANING

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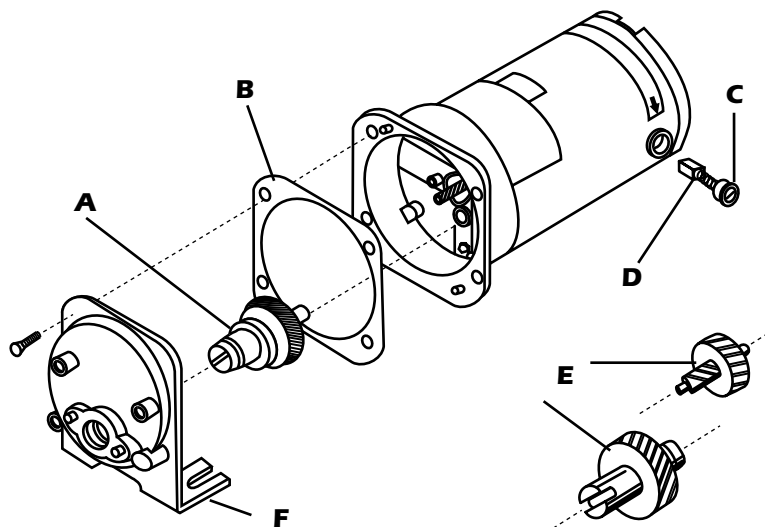
SPECIFICATIONS

Operating Temperature:	0° to 40°C (32° to 104°F)
Storage Temperature:	-45° to 65°C (-49° to 149°F)
Chemical Resistance:	ABS plastic case with a polyester label, gold, irridite-coated aluminum chassis and zinc-plated screws. All materials withstand standard cleaning solvents.
Line Voltage Limits:	90–130 V or 190–260 V, 50–60 Hz
Power Output:	0.075 kW (0.1 hp)
Maximum Current:	
115 V Units:	Controller: 2.2A, shorted output conditions
230 V Units:	1.1A, shorted output conditions
Start/Stop—Local Control:	
Input Voltage:	
Function Disable:	15V DC Typ.
Function Enable:	0.8V DC Max.
Input Current:	
Function Disable:	100 µA Max. leakage
Function Enable:	1.5mA Max.
Maximum Torque:	
300, 600 rpm Units:	13 kg•cm (180 oz-in)
100 rpm Units:	26 kg•cm (360 oz-in)
Speed Regulation:	
Line:	±1%
Load:	±2%
Speed Drift:	±10%
Enclosure Rating:	IP23 per IEC529
Humidity (non-condensing):	10% to 90%
Altitude:	Less than 2000 m
Display:	Green LED
Dimensions (L × W × H):	292 mm × 203 mm × 184 mm (11-1/2 in × 8 in × 7-1/4 in)
Weight:	5 kg (11 pounds)
Remote Control	
4–20 mA Input:	250 ohms typical input impedance referenced to signal ground. 4 mA, (Stop); 20 mA, (Full Speed) Accuracy: 3% F.S., Linear Resolution Overload Capability: 10V or 40 mA max.
2–10 V Input:	10 K ohms typical input impedance referenced to signal ground. 2 V, (Stop); 10 V, (Full Speed) Accuracy: 3% F.S., Linear Resolution Overload Capability: 15V max.
Compliance:	
115V	115V: UL508, CSA C22.2, No. 14-M91
230V	230V (For CE Mark): EN61010-1/A2: 1995 (EU Low Voltage Directive) and EN61326-1/A1: 1998 (EU EMC Directive)
Pollution Degree:	Pollution Degree 2 per IEC 664 (Indoor usage—lab, office)
Installation Category:	Installation Category II per IEC 664 (Local level—appliances, portable equipment, etc.)

REPLACEMENT PARTS AND ACCESSORIES; SPECIFICATIONS

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WARRANTY

Use only MASTERFLEX® precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is noted on the Warranty Card.

PRODUCT RETURN

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

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Vernon Hills, Illinois U.S.A. 60061-1844
Barrington, Illinois U.S.A. 60010-2392